



CARFIL QUALITY
YOUR SUPPLIER FOR ALL LABORATORY NEEDS

68U400 Electrode Angle Calibrator



This calibrator can correct the electrode position to any angle quickly and accurately in stereotaxic surgery. It can coordinate the position of the horizontal and vertical directions, the angle can be selected and move to the correct horizontal coordinate position, and the manipulator, which was placed in the dovetail arm, can move the tip of the electrode to the targeted coordinates in the vertical direction.

The horizontal slide has scales with 0,1mm accuracy and the zero point of it is correlated with that of the ear bars. When the horizontal slide is set at zero, the horizontal and vertical slide bar is fixed at zero position.

This calibrator is applicable only to the 68U000 series stereotaxes.



PAVAN SERVICE BVBA

Beyntellus 3 - B-2360 Oud-Turnhout - Tel. +32 (0)14 45 13 10 - admin@carfil.be - BTW BE 0420.681.872

www.carfil.be