




CARFIL QUALITY
YOUR SUPPLIER FOR ALL LABORATORY NEEDS

UNO Euthanasia Unit for IVC Cage

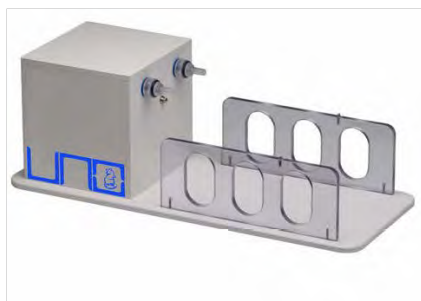

Features

A docking station, enabling human- and animal-welfare-compliant, intuitive and reliable euthanasia for mice in their home IVC. Just dock your IVC cage into the unit and the euthanasia process will start.

The UNO Euthanasia Unit for IVC-Cage has to be connected to an electrical wall outlet AND to a CO₂ cylinder or in-house CO₂ distribution system. The inlet pressure to the unit needs to be set at 0,5 bar via an external pressure regulator (not included).

When the IVC cage is correctly docked the unit ensures an even distribution of CO₂ and thus a homogeneously increasing concentration of CO₂ within the IVC cage. If the IVC cage is prematurely removed from the unit, the gas supply stops and the cycle ends prematurely.

The mice are euthanised in accordance with the AVMA Guidelines for the Euthanasia of Animals (2013), the German Animal Protection Act and Animal Welfare Regulation Governing Experimental Animals (2013).

The cycle begins when the CO₂ starts flowing into the unit and lasts 6 minutes to completion.

The UNO Euthanasia Unit for IVC-Cage is available for the following IVC cages:

- Allentown XJ
- Allentown Nexgen
- Tecniplast Blueline 1145
- Tecniplast Blueline 1284
- Tecniplast Blueline 1285 (L-type II NEXT)
- Tecniplast Greenline GM 50

Ordering number

180000201 + cage type

Euthanasia Unit IVC cage



PAVAN SERVICE BVBA

Beyntellus 3 - B-2360 Oud-Turnhout - Tel. +32 (0)14 45 13 10 - admin@carfil.be - BTW BE 0420.681.872

www.carfil.be