



CARFIL QUALITY
YOUR SUPPLIER FOR ALL LABORATORY NEEDS

Controlled Heating System Control Unit-MS/CU-MS



Control Unit-M/S

The **control Unit M/s** can like the Control Unit 02 be connected to a heating device and a temperature probe. However the interaction between the 2 connections allows you to **set the required animals rectal temperature**. The rectal probe measures the animals temperature and feed back is given to the controles of the heating device to keep the animals temperature as constant as possible. A safety feature is built-in to prevent over-heating in case the rectal probe is not in the animal. The heating element will not heat up over 42°C (also an acceptable norm for human skin).

With the Control Unit M/S you have a **self-regulating system** and do not need to manually adjust the temperature settings during the procedure.

Ordering number

Bestelnummer	
180000124	Control Unit MS
180000028	Flat Plate / 34,5 x 23 cm
180000128	Flat Plate / 18,5 x 11 cm
180000027	Half-Pipe Heating Unit
180000029	Tail Heating Unit
180000066	Temperature Probe - Mouse
180000026	Temperature Probe - Rat



PAVAN SERVICE BVBA

Beyntellus 3 - B-2360 Oud-Turnhout - Tel. +32 (0)14 45 13 10 - admin@carfil.be - BTW BE 0420.681.872

www.carfil.be