



CARFIL QUALITY
YOUR SUPPLIER FOR ALL LABORATORY NEEDS

Double 3- and 4-way valves

Eigenschappen



To be used for simultaneously changing the direction of both the **SUPPLY** of the anaesthetic gas mixture **AND** the **ECHAUST** of the anaesthetic gas mixture. The risk of making mistakes when the gas is supply and gas exhaust have to be changed individually from f.i. the Induction box to the face mask, is reduced substantially bij using these valves.

3-Way Valve

By turning the knob from the right to the left, both supply and exhaust are being directed from the face mask to/from the Induction box.

4- Way Valve

- A.. By turning the knob to the left, both supply and exhaust are being directed to/from the induction box.
- B.. By turning the knob to the center, both supply and exhaust are being directed to/form the intubation aid.
- C.. By turning the knob to the right, both supply and exhaust are being directed to/from the UMV-03

Bestelnummer

180000159	Double 3-way valve
180000259	Double 4-way valve



PAVAN SERVICE BVBA

Beyntellus 3 - B-2360 Oud-Turnhout - Tel. +32 (0)14 45 13 10 - admin@carfil.be - BTW BE 0420.681.872

www.carfil.be